# Knick >

# The requirements for lab measurements become stricter every day. Quality assurance and measurement documentation in accordance with GLP are a must in many areas.

With its numerous safety functions and record printouts at keystroke, the 703 Laboratory Conductivity Meter considerably simplifies this work for you.

#### **Fullcheck**

automatically checks the device functions during power-on. Also during operation, a complete instrument check can be carried out at a single keystroke. Here, also display and keypad are checked besides the electrical characteristics.

#### **Record printouts**

With record printouts of the device self-test, the calibration, and the parameter settings, it is possible (as part of quality management to ISO 9000 and GLP) to document the operability and the regular maintenance and calibration of the meter.

#### Sensoface

Sensoface monitors the sensor and measuring equipment and provides information on sensor selection and handling. It reports clock memory loss and requests regular checks in accordance with GLP.

## **Calibration**

Unknown cell constants can easily be determined with a standard calibration. The meter automatically takes the TC of the calibration solution into consideration, calculates the cell constant and displays it. Of course, a known cell constant can also be entered

# 703 Laboratory Conductivity Meter



directly.

#### **Analog output**

The galvanic isolation of the recorder output prevents the measured values from being influenced by the connected peripherals. Measurement continues unimpaired.

#### **EMC**

EMC design protects the meter from electromagnetic interferences, ensuring reliable measurement results even under unfavorable conditions.

This makes the Model 703 the first laboratory conductivity meter that completely fulfills the EMC recommendations of NAMUR.

The Model 703 offers a wide range of practical features to meet the numerous requirements of everyday measuring tasks.

# Automatic switchover to 4-electrode or 2-electrode operation

With the Model 703 you can use either 4-electrode or 2-electrode sensors. The measuring input is automatically switched to the appropriate operating mode.

# Temperature compensation manual or automatic

Temperature compensation takes place either automatically with Pt 1000/NTC 30 kOhm temperature probes or manually.

#### Standard RS 232 interface

Via the standard RS 232 interface your data can be immediately processed by a computer. Even direct output to a printer is no problem.

### GLP records at the press of a key

Records of the parameter setting, calibration, and device diagnostics can be output directly to a printer. This provides you with comprehensive GLP-compatible documentation at the press of a key.

# Automatic adjustment of display range

The meter automatically selects the display range with the greatest possible resolution. Of course, the desired display range can also be specified manually.

# Easy-to read LED display for two measured values

The large, bright LED display allows simultaneous readout of two measured values, such as conductivity and temperature. The 14-segment display can show alphanumeric characters.

## Double insulation provides electrical safety in wet locations

The well-designed enclosure has proved successful in practical use. A waterproof membrane keyboard and drain grooves protect the meter from moisture. The robust, stainless steel covered enclosure resists even strong mechanical stress.

# **Conductivity Measurement**







Knick >



Measurement ranges from  $0.000 \mu S/cm$  to 2000 mS/cm

<1.000  $\mu$ S/cm ... >1000 mS/cm with one sensor

Records for QM documentation to ISO 9000 and GLP

Calibrated analog recorder output, galvanically isolated

Sensoface monitoring of sensor and measuring equipment

Automatic calibration with standard solutions

EMC to NAMUR

RS 232 interface for computer and printer



Two measured value displays, simultaneous

Self-contained clock

Liquid-proof membrane keypad

Robust enclosure

IP 54 protection

3-year warranty



# Warranty vears.

Defects occurring within 3 years from delivery date shall be remedied free of charge at our plant (carriage and insurance paid by sender). Sensors and accessories: 1 year




# **703 Laboratory Conductivity Meter**

## Keypad

Exit function and return to measuring mode

Print currently measured values or function data Select line, edit value or select variable Select parameter or position

Select line, edit value or select variable

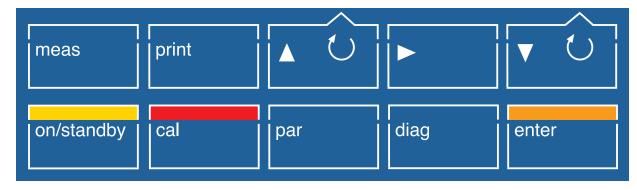

On/off (standby)

Activate calibration

Activate parameter setting

Activate diagnostics

Take over value or entry

## **■** Record printouts

Records of parameter setting, calibration and diagnostics are particularly helpful for QM documentation to ISO 9000 and GLP.

The records can be printed out directly to any commercially available printer with serial port.

Serial Number: Software Version: Hardware Version: Options:	01100329 1.3 01 No
Last Calibration:	19.03.03 10:03
Data Entry	OF DESCRIPTION OF SECURE AS A SECURE OF SECURE AS A
Cell System Data Cell Constant:	1.240/cm

Knick 703 Diagnost	tics 19.03.0
Serial Number: Software Version: Hardware Version: Options:	01108329 1.3 01 No
Last Fullcheck:	19.03.03 09:55
RAM: PROM: EEPROM: Output: Amplifier: Accumulator: Display: Keys:	-ok- -ok- -ok- -ok- -ok- - tested- -ok-
Sensoface(++/oo/)	
Cell Range: TC Temperature GLP Timer: Accumulator: Date Check:	÷ ÷ ÷ ÷ ÷ ÷ ÷ ÷ + ÷ + ÷ + † † † † † † †
nate rueck:	***

Serial Number: Software Version: Hardware Version:	01108329 1.3 01
Options : 	No
Manual Temperature:	25.0CEL
Temperature Compensat	ion: On
Temperature Coefficie	nt: 2.10%/K
Reference Temperature	
Sensoface:	<u>On</u>
Range:	Fixed
Cal-Solution:	MaCl 0.1Mol
GLP Timer:	0h
Recorder Output: Baud Rate:	1 mU/mS 48200
Data Bits/Parity:	7 Even
Protocol:	No Everi
Interface:	Printer
Printer Timer:	0.0min
Time:	10:47
Date:	19.03.
vace: Year:	17.05. 2003
1.001 .	7609

# Knick >


#### ■ Fullcheck device self-test

For the self test the sensor is automatically switched off and the input switched over to a reference resistor. The conductivity measuring circuit is automatically checked internally.

**Conductivity Measurement** 

The microcontroller sends defined voltage steps to the recorder output. These are measured with the A/D converter and compared with a highly accurate reference voltage.

This means, a complete test of the signal path is implemented with a conductivity meter for the first time. In addition, all memories, the display, and the keypad are tested.

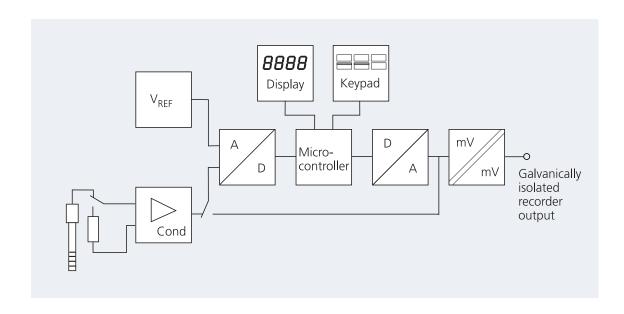

**Conductivity Measurement** 

# **703 Laboratory Conductivity Meter**

# **■** Specifications

Equipment	Meter with power cord, without sensor		
Ranges	Cond: 0.000 9.999 μS/cm; 00.00 99.99 μS/cm; 000.0 999.9 μS/cm; 0.000 99.99 mS/cm; 00.00 99.99 mS/cm; 0000 999.9 mS/cm; 0000 2000 mS/cm; auto-ranging or manual preset*);  °C: –50.0 +150.0		
Display	Alphanumer. 2 x 4-digit, 14-segment LED, character height 13 mm, measurement symbols: 20 °C, 25 °C, μS/cm, mS/cm, %/K, °C, Time, 3 Sensoface icons inform on sensor and measuring equipment (GLP)		
Measuring cycle	Approx. 1.5/sec		
Measuring frequencies	Approx. 40 Hz to 2 kHz, automatic adjustment by conductance		
Resolution	Γ		
Accuracy <sup>1)</sup>	Cond: <0.5 % meas. value. ±2 counts °C: 0.3 K		
Reproducibility <sup>1)</sup>	< 0.1 % meas. val.		
Temperature compensation	−50 +150 °C; Pt 1000/NTC 30 kOhm (autom. selection) or manual, Linear TC characteristic 0.00 +9.99 %/K, ref. temperature 20 °C/25 °C selectable		
Adm. cell constant	0.001 199.9 cm <sup>-1</sup> , adjustable		
Sensor standardization	Operating modes – Automatic by determining the cell constant with NaCl or KCl solution Calibration solutions: KCl 0.01 mol/l; 0.1 mol/l; 1 mol/l; NaCl 0.01 mol/l; 0.1 mol/l; saturated – Direct entry of cell constant		
Monitoring of sensor and equipment (GLP)	Sensoface provides information:  - for selection of 2-electrode sensors  - on too great a difference between reference and measuring temperature  - for handling of 4-electrode sensors  - on clock memory loss  - in case of irregular checking of measuring equipment  Optical display: good/average/poor		
Device self-test	Test of measuring electronics including recorder output, segment and keypad test, RAM, EPROM, and EEPROM test during diagnostics, automatic short check at power-on		
GLP records (ISO 9000)	Parameter settings, calibration, device diagnostics		
Recorder output	Galvanically isolated (isolation voltage: 40 V DC, 20 V AC) Cond: 1 mV/ $\mu$ S · cm $^{-1}$ ; 1 mV/mS · cm $^{-1}$ ; °C: 10 mV/°C user-defined for printer control		
nterface	RS 232 without control lines, galvanically isolated (isolation voltage: 40 V DC, 20 V AC), user-definable as printer or computer interface,  Baud rate: 600 / 1200 / 2400 / 4800 / 9600*) Data bits/parity: 7/Even, 7/Odd, 8/No*)  Protocol: none, XON/XOFF*) Stop bits: 1		
Software	Control of the Model 703 Laboratory Conductivity Meter is integrated in the automation software for lab meters "labworldsoft" (Fisher Scientific) for display and control of device		

functions for Version 4.0 or higher.

**Conductivity Measurement** 

Clock 	Real-time clock with date, self-contained
Calibration data storage	Automatic storage of cell constant and calibration procedure with time and date stamp, self-contained
Oata retention	Parameters, statistics, and factory settings: >10 years (EEPROM)  Clock: reserve power >1 year (battery-backed)
rotection against electrical hock	Protective separation as defined in DIN 57100 / VDE 0100 Part 410 and DIN VDE 0106 Part 101, power supply against all other inputs and outputs, in accordance with the NAMUR recommendation "Extra-low voltage circuits with protective separation"
MC directive	89/336/EEC
tandards	EN 61326 / VDE 0843 Part 20: 2002-3
Ambient temperature Storage and transport temp	0 +45 °C -20 +70 °C
Power supply	230 V –15 % +10 %, 48 62 Hz, <10 VA, Option 363: 115 V AC
ensor connection	The meter allows connection of any 2-electrode sensors with banana plug.  Special diode plug for 4-electrode sensors.
Enclosure	Glass-reinforced polyamide 12, stainless steel cover, IP 54 protection, prepared for connecting ZU 6954 attachable stand
Dimensions (W x H x D)	244 x 95 x 255 mm
Veight	Approx. 2 kg

Conductivity Measurement

## **■** Specifications Accessories

Printer	Order No.: ZU 0244		
Туре	Matrix printer		
Interface	Serial RS 232 port		
Paper	Standard paper, width: 57.5 mm (2.25 inches)		
Baud rate	4800 bauds		
Data bits	7, 1 stop bit		
Parity	Even		
Protocol	No No		
Power supply	230 V AC ±10 %		
Dimensions (W x H x D)	197 x 73 x 153 mm		
Weight	Approx. 1.2 kg (incl. power pack)		
Stand	Order No.: ZU 6954		
Material	Pillar: anodized aluminum; carriage and base: polyamide 12 glass reinforced; Beaker stop, vertical stop, and electrode clasp: stainless steel		
Carriage stroke	190 mm		
Clamping possibilities Stop for sample beakers	2 x 12 ±0.5 mm; 1 x 4 14 mm; 1 x 6 16 mm from Ø 30 150 mm		
Beaker height	Г — Г — Г — Г — Г — Г — Г — Г — Г — Г —		
Dimensions (W x H x D)	130 x 300 x 145 mm		
Weight	Approx. 410 g		
Plug-in power pack for immersion stirrer	Order No.: ZU 6956		
Power supply	230 V AC –15 % +6 % <8 VA		
Cable length	2 m		
Weight	Г Арргох. 380 g		
Immersion stirrer	Order No.: ZU 6955		
Г————————————————————————————————————	Enclosure: PVC; impeller and shaft: stainless steel		
Dimensions	Unit: 250 x Ø 25/12 mm; impeller: Ø 12 mm; immersion depth: approx. 90 mm		
Weight	Г		

#### Conductivity sensors for lab and portable meters

**Conductivity Measurement** 

SE 202 2-electrode sensor with integrated temperature probe (NTC 30 kOhm) and flow cell. For measurement in low-conductivity solutions such as ultrapure water and boiler feed water, e.g. for monitoring water desalination plants.

SE 204 4-electrode sensor with integrated temperature probe (NTC 30 kOhm). For measurement in natural waters such as surface water or drinking water, in aqueous solutions such as acid and alkaline solutions and for salinity determination of sea water.

With the ZU 6985 4-electrode sensor from Knick, a lab-quality universal conductivity sensor is available. The sensor operates reliably over a broad range from <1.00  $\mu\text{S/cm}$  to >1000 mS/cm. It is equipped with a quick-reacting Pt 1000 temperature probe. It is provided with a glass/platinum measuring system with an easy-to-replace KPG tube. It is simple to clean and requires no platinization.

#### **■** Specifications Conductivity sensors

Conductivity sensors	SE 202	SE 204	ZU 6985
Number of electrodes	2	4	4
Body Electrode material	Stainless steel 1.4571 Stainless steel 1.4571	Epoxy, black Graphite	Glass Platinum, bare
Body length	120 mm	120 mm	110 mm
Body diameter	12 mm	15.3 mm	Tube 16 mm
Temperature probe	NTC (30 kOhm): -5 +100 °C	NTC (30 kOhm): -5 +100 °C	Pt 1000: −20 +100 °C
Immersion depth	Min.: 30 mm, max.: total length incl. Cable	Min.: 36 mm, max.: total length incl. Cable	Min. 60 mm Max. 80 mm
Pressure resistance	2 bars	2 bars	2 hars
Cell constant	$0.100 \text{ cm}^{-1} \pm 2 \%$	0.475 cm <sup>-1</sup> ±1.5 %	1.19 cm <sup>-1</sup> ±1 %
Ranges	0.01 200 μS/cm	1 μS/cm 500 mS/cm	1 μS/cm 1000 mS/cm
Remarks	Incl. flow cell	-	
Order No.	SE 202	SE 204	ZU 6985
The state of the s			

# ■ Product line Laboratory conductivity meters and conductivity sensors

Lab Conductivity Meter	703	Order No.
Options	Unit with power cord, without sensor	703
Power supply	115 V AC	363
	F	
2-electrode sensor	With stainless steel body incl. flow cell (ZU 0298 adapter required)	SE 202
4-electrode sensor	With epoxy body (ZU 0298 adapter required)	SE 204
4-electrode sensor	With glass body	ZU 6985

Order No.

## ■ Product line Calibration solutions, spare parts, and further accessories

**Conductivity Measurement** 

190	Temperature probe	Pt 1000	ZU 6959
The state of the s	Conductivity standard	For determination and checking of cell constants.  1 ampoule for producing 1000 ml 0.1 mol/l NaCl solution (12,88 mS/cm)	ZU 6945
		For determination and checking of cell constants. Conductivity 12.88 mS/cm $\pm 1$ % (0.1 mol/l KCl), 250 ml solution, ready for use	ZU 0348
		For determination and checking of cell constants. Conductivity 1413 $\mu$ S/cm $\pm$ 1 % (0.01 mol/l KCl), 250 ml solution, ready for use	ZU 0349
		For determination and checking of cell constants. Conductivity 147 $\mu$ S/cm $\pm$ 1 %, 500 ml solution, ready for use	ZU 0702
		For determination and checking of cell constants. Low conductivity 15 $\mu$ S/cm $\pm$ 5 %, 300 ml solution, ready for use	ZU 0350
0	KPG tube	For ZU 6985 4-electrode sensor, incl. O-ring	ZU 0180
	Replacement flow cell	For SE 202 2-electrode sensor	ZU 0284
	Adapter	For connecting the SE 202 and SE 204 sensors to the 703 Laboratory Conductivity Meter	ZU 0298



Attachable stand

Besides the immersion stirrer, the attachable stand can hold three sensors of any kind. The adjustable stops prevent damage of sensor and beaker glass. Timeconsuming adjustment during sample changes has been eliminated. An integrated cable duct does away with the "spaghetti cables" on your benchtop. For ZU 6955 immersion stirrer and three sensors, directly connected to the meter.

ZU 6954

continued – **Product line** Calibration solutions, spare parts, and further accessories



Immersion stirrer

The immersion stirrer reduces sensor response time for measurement and calibration. Precision measurements to DIN 19268 even require stirring. To prevent splattering of test liquid, the stirrer automatically stops as the carriage moves up. The stirrer is supplied via the ZU 6956 plug-in power pack.

Order No.

ZU 6955

92



## **■** Product line Accessories

			Order No.
	Plug-in power pack	For immersion stirrer	ZU 6956
	Interface cable	For meter – computer connection (special EMC cable)	ZU 0152
1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1	Lab printer	With the Lab Printer, you can document your measured values either at the press of a key or timer-controlled. Also records for QM documentation to ISO 9000 and GLP can be printed out with a single keystroke. The printer is equipped with a replaceable ribbon cartridge and prints on standard paper. It is connected to the 765 Laboratory pH Meter or the 703 Laboratory Conductivity Meter via interface cable.	ZU 0244
	Interface cable	For meter – printer connection	ZU 0245
I	Printer paper	For ZU 0244 Lab Printer, 5 rolls	ZU 0249
The state of the s	Ink ribbon	For ZU 0244 Lab Printer, 5 ribbons	ZU 0250