# Itton/L

Compact, in-compartment microscope for the FT/IR-4000/6000 Series




U/S.CO



measurements of samples approaching 20 microns.

The Irtronµ sample compartment microscopy system is designed to The Irtronµ can be operated with the integrated touch panel or via a PC provide affordable analysis of microscopic samples with the high using the Spectra Manager™ II software interface. The sample is performance features of an external FT-IR microscope accessory. The observed with the integrated color CCD camera and 5 inch TFT color Irtronu offers unprecedented convenience and ease of use, compatible LCD monitor. JASCO's unique ATOS (Aperture Through Optical with the JASCO FT/IR-4000 and 6000 Series FT-IR instruments. The System) provides simultaneous viewing of the sample image and microscope accessory installs into the spectrometer sample sampling location specified by the aperture. The ATOS combined with compartment in seconds without optical alignment. The state-of-the-art the SmartMonitor™ feature allows observation of the sample while optical design guarantees high throughput for highly sensitive collecting spectra in real-time using the monitor mode. An option within the Spectra Manager control package allows auto-saving of the collected spectra as well as the sample image(s).

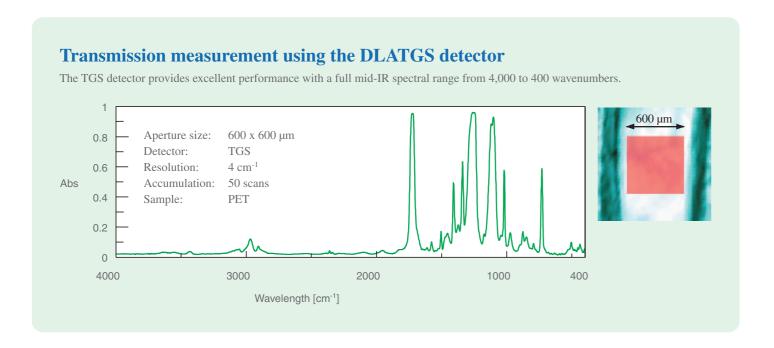


- Transmittance, reflectance and ATR (option) measurement modes
- **DLATGS, MCT or NIR optimized detector**
- Unique SmartMonitor™ function allows sample observation during spectrum preview
- Five inch TFT LCD monitor for sample observation
- Optional sampling accessories for liquid and powder samples
- Affordable, easy-to-use microscopy system



#### Transmission measurement with the MCT detector

The MCT detector provides excellent sensitivity and sample measurements approaching 20 microns.



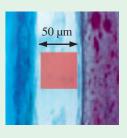






## **SmartPurge** TM

An optional purge casing can be used to eliminate interference from CO<sub>2</sub> and water vapor bands.

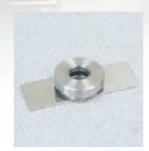

# Liquid sample holder

The liquid sample holder includes three standard spacers; 0.1, 0.05 and 0.025 mm pathlengths for liquid sample analyses.



## **ATR** objective

Three types of optional ATR objectives are available, selecting from ZnSe, Ge, and Diamond crystal elements.



# Powder sample holder

The powder sample holder can be used for diffuse reflectance measurements of powdered samples mixed with KBr. Up to five samples can be loaded onto the holder.



### **Pressure sensor**

An optional pressure sensor (recommended) to protect the ATR objective is also available.



## **SliceMaster** TM

SliceMaster TM is a handy compact instrument that can easily create thin sections by cutting film-type samples. The SliceMaster is a low-cost cutting tool with capabilities comparable to a general purpose microtome.



• Specifications are subject to change without notice.

## JASCO INTERNATIONAL CO., LTD.

4-21, Sennin-cho 2-chome, Hachioji, Tokyo 193-0835, Jápan
Tel: +81-426-66-1322 Fax: +81-426-65-6512 Internet: http://www.jascoint.co.jp/english/index.html
Australia, China, Hong Kong, India, Indonesia, Iran, Korea, Malaysia, New Zealand,
Pakistan, Philippines, Russia, Singapore, South Africa, Taiwan, Thailand

#### JASCO INCORPORATED

8649 Commerce Drive, Easton, Maryland 21601-9903, U.S.A
Tel:+1-800-333-5272 Tel:+1-410-822-1220 Fax:+1-410-822-7526 Internet:http://www.jascoinc.com
Canada, Costa Rica, Mexico, Puerto Rico, Argentina, Brazil, Chile, Colombia, Paraguay, Peru, Uruguay

#### JASCO EUROPE S.R.L.

Via Cadorna 1, 23894 Cremella (Lc), Italy

Tel: +39-039-9215811 Fax: +39-039-9215835 www.jasco-europe.com

 $\textbf{JASCO Deutschland} \ www.jasco.de, \ \textbf{JASCO UK} \ www.jasco.co.uk, \ \textbf{JASCO France} \ www.jascofrance.fr,$ 

JASCO Benelux www.jasco.nl, JASCO Spain www.jasco-spain.com,

Austria, Finland, Greece, Hungary, Poland, Portugal, Romania, Switzerland, Algeria, Cyprus, Egypt, Israel, Jordan, Kuwait, Lebanon, Morocco, Saudi Arabia, Syria, Tunisia, Turkey, U.A.E.

