


# Extraction of 45 Multi Class Drugs of Abuse from Hydrolyzed Urine Using ISOLUTE® SLE+ prior to LC-MS/MS Analysis

This application note describes a Supported Liquid Extraction (SLE) protocol for the extraction of various drugs of abuse from hydrolyzed urine prior to LC-MS/MS analysis.

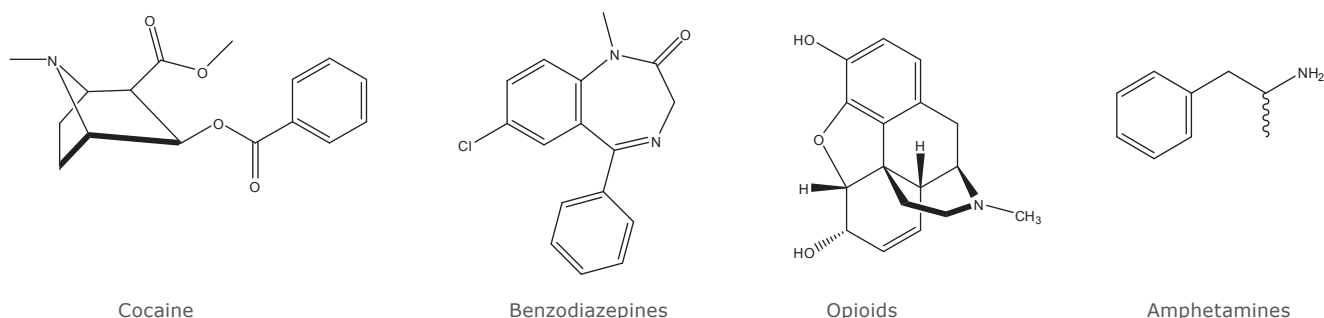

**Figure 1.** Example structures by class

## Introduction

The method described in this application note achieves high reproducible recoveries for a wide range of drugs of abuse and metabolites from hydrolyzed urine. Hydrolyzed urine was extracted using the ISOLUTE SLE+ 200 supported liquid extraction plate and 1 mL column formats using 100 and 500  $\mu$ L sample volumes, respectively. Sub ng/mL levels were achieved for all analytes using 100  $\mu$ L of urine on the fixed well plate format.

ISOLUTE® SLE+ products provide clean, rapid, robust, efficient, high throughput and automatable extraction solutions for a wide range of analytes.

## Analytes

Amphetamine, Methamphetamine, MDA, MDMA, MDEA, Methadone, EDDP, Mephedrone, Morphine, Hydromorphone, Oxymorphone, Dihydrocodeine, Oxycodone, Hydrocodone, Codeine, 6-MAM, Cocaine, Benzoylecgonine, Fentanyl, Norfentanyl, Ketamine, Norketamine, Buprenorphine, Norbuprenorphine, 7-amino-flunitrazepam, 7-amino-clonazepam, Nitrazepam, Flunitrazepam, Clonazepam,  $\alpha$ -OH-alprazolam,  $\alpha$ -OH-triazolam, Oxazepam, Estazolam, Temazepam, Alprazolam, Lorazepam, 2-OH-ethyl-flurazepam, Triazolam, Nordiazepam, Diazepam, Midazolam, Flurazepam, Zaleplone, Zopiclone, Zolpidem.

## Sample Preparation Procedure

|                         |   |
|-------------------------|---|
| <b>SLE+ Format</b>      | ISOLUTE® SLE+ 200 supported liquid extraction plate, part number 820-0200-P01 or ISOLUTE SLE+ 1 mL sample volume supported liquid extraction columns, part number 820-0140-C.   |
| <b>Urine Hydrolysis</b> | <p>Take 1 mL urine and spike with internal standard (10 µL of Amphetamine-d3, Morphine-d3, Diazepam-d5, parent concentration 250 ng/mL; BZE-d3 and 6-MAM-d3, metabolite concentration 25 ng/mL). Add 1 mL of 100 mM ammonium acetate buffer (pH5) and 50 µL β-Glucuronidase enzyme (equivalent to approximately 4500 U/mL of urine).</p> <p>Hydrolyze at 60 °C for 2 hours.</p> |
| <b>Post Hydrolysis</b>  | <p>Cool the hydrolyzed urine and add 10 µL of concentrated ammonium hydroxide (28–30% stock, aq) per mL of urine.</p> <p>Vortex mix thoroughly.</p>   |

## Supported Liquid Extraction

|                       |  |
|-----------------------|--|
| <b>Sample Loading</b> | <p><b>ISOLUTE SLE+ 200 µL plate:</b> Load pre-treated sample (200 µL) to the 96 well-plate followed by a pulse of vacuum or positive pressure to initiate flow. Leave to absorb for 5 minutes.</p> <p><b>ISOLUTE SLE+ 1 mL columns:</b> Load pre-treated sample (1 mL) to the columns followed by a pulse of vacuum or positive pressure to initiate flow. Leave to absorb for 5 minutes.</p>  |
| <b>Elution</b>        | <p><b>ISOLUTE SLE+ 200 µL plate:</b> Apply 1 mL of dichloromethane/isopropanol (95/5, v/v) and allow to flow under gravity for 5 minutes. Pull through the remaining solvent with vacuum or positive pressure for 10–20 seconds.</p> <p><b>ISOLUTE SLE+ 1 mL columns:</b> Apply 2.5 mL of dichloromethane/isopropanol (95/5, v/v) and allow to flow under gravity for 5 minutes.</p> <p>Apply a second 2.5 mL aliquot of dichloromethane/isopropanol (95/5, v/v) and allow to flow under gravity for 5 minutes. Pull through the remaining solvent with vacuum or positive pressure for 10–20 seconds.</p> <p>Note: The addition of 100 µL of 50 mM HCl in methanol into the collection plate or culture tube is required prior to or post elution to stabilize multiple analytes (amphetamines, bath salts and ketamine) due to volatility issues during evaporation.</p> |
| <b>Post Elution</b>   | Evaporate to dryness at 40 °C in a stream of air or nitrogen using a SPE Dry.  |
| <b>Reconstitution</b> | <p><b>ISOLUTE SLE+ 200 µL plate:</b> Reconstitute in 200 µL of reconstitution buffer. Cap with a sealing mat and vortex gently.</p> <p><b>ISOLUTE SLE+ 1 mL columns:</b> Reconstitute in 200 µL of reconstitution buffer and vortex gently.</p>  |

## UPLC Conditions

|                     |   |
|---------------------|---|
| <b>Instrument</b>   | Waters ACQUITY UPLC 20 $\mu$ L Loop                                 |
| <b>Column</b>       | ACQUITY UPLC BEH C18 column (1.7 $\mu$ m, 100 x 2.1 mm id)          |
| <b>Mobile Phase</b> | A: 2 mM ammonium acetate (aq)<br>B: 2 mM ammonium acetate /methanol |
| <b>Flow Rate</b>    | 0.3 mL/min  |

**Table 1.** Gradient Conditions

| Time | % A | % B | Curve |
|------|-----|-----|-------|
| 0    | 90  | 10  | 1     |
| 10   | 10  | 90  | 6     |
| 11.4 | 10  | 90  | 6     |
| 13.4 | 90  | 10  | 1     |

**Curve 1.** Conditions in line initiated immediately once previous time passed. i.e. 90:10 resumed at 11.4 minutes.

**Curve 6.** Linear Gradient

|                           |   |
|---------------------------|---|
| <b>Injection Volume</b>   | 10 $\mu$ L (partial loop with overfill) |
| <b>Sample Temperature</b> | 20 °C                                   |
| <b>Column Temperature</b> | 40 °C                                   |

## Mass Spectrometry Conditions

|                                |  |
|--------------------------------|--|
| <b>Instrument</b>              | Waters Premier XE triple quadrupole mass spectrometer equipped with an electrospray interface for mass analysis. |
| <b>Desolvation Temperature</b> | 450 °C   |
| <b>Ion Source Temperature</b>  | 120 °C   |

Positive ions acquired in the multiple reaction monitoring (MRM) mode:


**Table 2.** MRM Conditions

| Compound                | MRM Transition | Cone Voltage (V) | Collision Energy (eV) |
|-------------------------|----------------|------------------|-----------------------|
| Amphetamine             | 136.0 > 118.9  | 16               | 9                     |
| Amphetamine-d5          | 141.0 > 123.9  | 16               | 9                     |
| Methamphetamine         | 150.0 > 90.9   | 22               | 17                    |
| MDA                     | 180.1 > 105.0  | 16               | 23                    |
| MDMA                    | 194.1 > 163.0  | 20               | 13                    |
| MDEA                    | 208.2 > 163.0  | 22               | 13                    |
| Hydromorphone           | 286.2 > 185.1  | 44               | 29                    |
| Morphine                | 286.2 > 201.0  | 42               | 25                    |
| Morphine-d3             | 289.2 > 201.0  | 42               | 25                    |
| BZE                     | 290.1 > 168.0  | 30               | 18                    |
| BZE-d3                  | 293.1 > 171.0  | 30               | 18                    |
| Oxymorphone             | 302.2 > 198.1  | 34               | 37                    |
| Dihydrocodeine          | 302.2 > 199.1  | 42               | 33                    |
| Oxycodone               | 316.2 > 241.2  | 34               | 27                    |
| Mephedrone              | 178.1 > 160.0  | 35               | 12                    |
| Norfentanyl             | 233.1 > 84.0   | 25               | 19                    |
| 7-amino-flunitrazepam   | 284.2 > 135.0  | 40               | 27                    |
| 7-amino-clonazepam      | 286.2 > 121.0  | 40               | 30                    |
| Hydrocodone             | 300.2 > 199.1  | 46               | 33                    |
| Codeine                 | 300.3 > 215.1  | 42               | 25                    |
| 6-MAM                   | 328.2 > 165.1  | 44               | 33                    |
| 6-MAM-d3                | 331.2 > 165.1  | 44               | 33                    |
| Cocaine                 | 304.2 > 182.0  | 30               | 20                    |
| Norketamine             | 224.1 > 124.9  | 20               | 23                    |
| EDDP                    | 278.2 > 234.2  | 26               | 30                    |
| Zaleplone               | 306.2 > 264.2  | 40               | 22                    |
| Zopiclone               | 389.2 > 245.1  | 20               | 17                    |
| Norbuprenorphine        | 414.3 > 101.0  | 55               | 42                    |
| Ketamine                | 238.1 > 124.9  | 25               | 27                    |
| Nitrazepam              | 282.2 > 236.1  | 40               | 25                    |
| Flunitrazepam           | 314.2 > 268.2  | 40               | 25                    |
| Clonazepam              | 316.1 > 270.1  | 40               | 25                    |
| $\alpha$ -OH-triazolam  | 359.1 > 331.1  | 45               | 26                    |
| Oxazepam                | 287.2 > 241.0  | 30               | 21                    |
| Estazolam               | 295.2 > 267.2  | 40               | 24                    |
| Temazepam               | 301.1 > 255.1  | 30               | 22                    |
| Zolpidem                | 308.2 > 235.1  | 45               | 35                    |
| Alprazolam              | 309.2 > 281.2  | 40               | 26                    |
| Methadone               | 310.2 > 265.2  | 26               | 15                    |
| Lorazepam               | 321.1 > 275.1  | 30               | 22                    |
| $\alpha$ -OH-alprazolam | 325.2 > 297.1  | 40               | 25                    |
| 2-OH-ethyl-flurazepam   | 333.2 > 109.0  | 40               | 27                    |
| Triazolam               | 343.0 > 308.1  | 45               | 27                    |
| Nordiazepam             | 271.1 > 139.9  | 40               | 28                    |
| Diazepam                | 285.2 > 154.0  | 40               | 27                    |
| Diazepam-d5             | 290.2 > 154.0  | 40               | 27                    |
| Midazolam               | 326.2 > 291.2  | 45               | 29                    |
| Fentanyl                | 337.3 > 105.0  | 35               | 40                    |
| Flurazepam              | 388.2 > 315.1  | 35               | 23                    |
| Buprenorphine           | 468.3 > 101.0  | 55               | 42                    |


## Results

### Recovery


Urine spiked with 2 ng of analytes (n=7), equating to 20 ng/mL or 4 ng/mL when extracting 100 or 500  $\mu$ L of urine, respectively. RSD's (n=7) ranged from 0.1–9.6%.

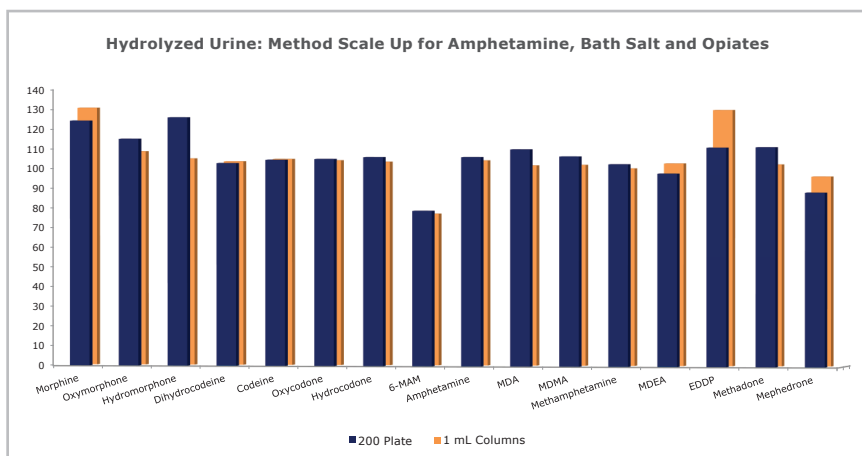

**Figure 2.** Recovery profile for amphetamines, bath salt and opiates from hydrolyzed urine using ISOLUTE® SLE+ 200 fixed well plates and 1 mL columns.

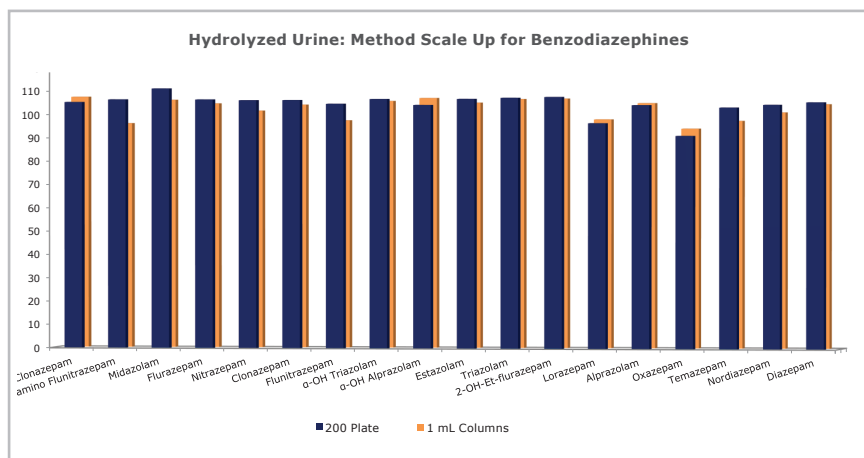


**Figure 3.** Recovery profile for benzodiazepines from hydrolyzed urine using ISOLUTE® SLE+ 200 fixed well plates and 1 mL columns.

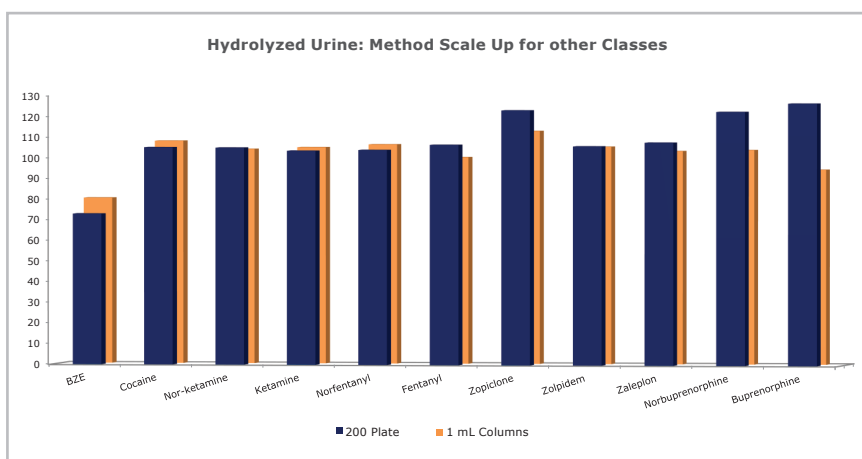


**Figure 4.** Recovery profile for multi class analytes from hydrolyzed urine using ISOLUTE® SLE+ 200 fixed well plates and 1 mL columns.

## Calibration Curves


200  $\mu$ L fixed well plate and 1 mL column processing: Calibration curves were generated using urine spiked at concentrations from 1–500 ng/mL. Good coefficients of determination were obtained for all analytes ( $r^2 > 0.99$ ). Quadratic function was observed at the top end of the calibration curve for many analytes. Dilution of these samples was performed to improve linearity.

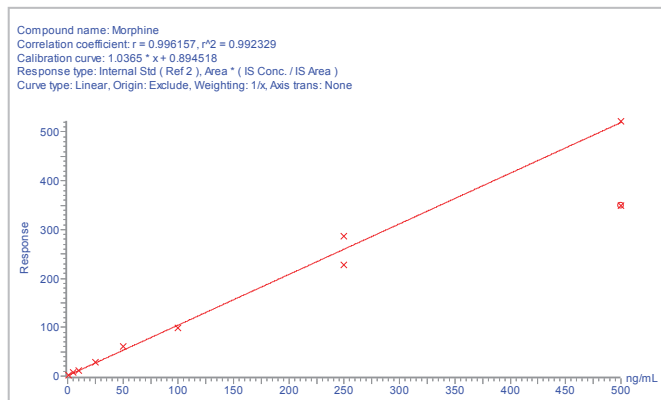


**Figure 5.** Calibration Curve for morphine using ISOLUTE® SLE+ 1 mL columns.

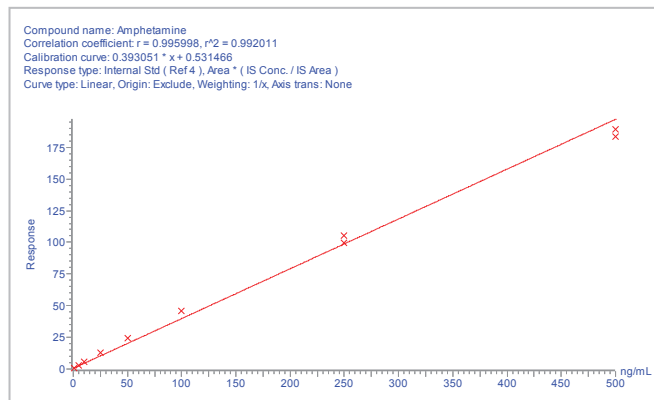

**Figure 6.** Calibration Curve for amphetamine using ISOLUTE SLE+ 1 mL columns.

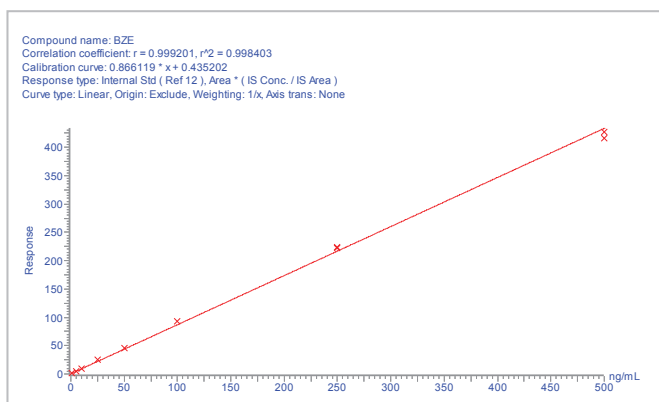

**Figure 7.** Calibration Curve for benzoylcegonine (BZE) using ISOLUTE® SLE+ 1 mL columns.

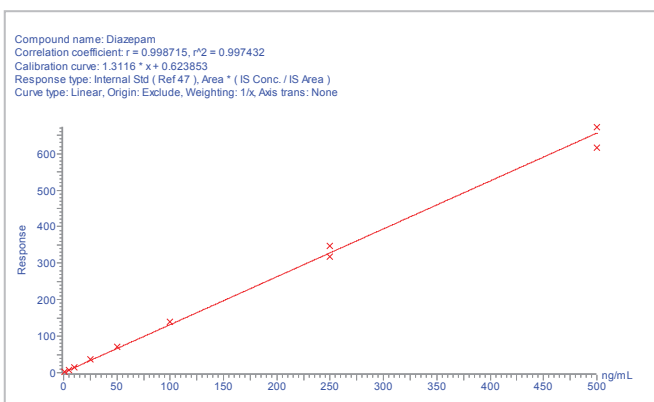

**Figure 8.** Calibration Curve for diazepam using ISOLUTE SLE+ 1 mL columns.

**Table 3.** Estimated LOQs based on S/N ratios from 1 ng/mL extracted samples are:

| Analyte               | Estimated LOQ<br>200 µL Plate ng/mL | Estimated LOQ 1 mL<br>Columns ng/mL |
|-----------------------|-------------------------------------|-------------------------------------|
| Amphetamine           | < 0.3                               | < 0.2                               |
| Methamphetamine       | < 0.15                              | < 0.1                               |
| MDA                   | < 0.2                               | < 0.1                               |
| MDMA                  | < 0.1                               | < 0.05                              |
| MDEA                  | < 0.2                               | < 0.1                               |
| Hydromorphone         | < 0.2                               | < 0.2                               |
| Morphine              | < 0.5                               | < 0.3                               |
| BZE                   | < 0.1                               | < 0.05                              |
| Oxymorphone           | < 0.5                               | < 0.5                               |
| Dihydrocodeine        | < 0.1                               | < 0.1                               |
| Oxycodone             | < 0.5                               | < 0.5                               |
| Mephedrone            | < 1                                 | < 0.75                              |
| Norfentanyl           | < 0.1                               | < 0.05                              |
| 7-amino-flunitrazepam | < 0.5                               | < 0.25                              |
| 7-amino-clonazepam    | < 0.5                               | < 0.25                              |
| Hydrocodone           | < 0.25                              | < 0.2                               |
| Codeine               | < 0.25                              | < 0.2                               |
| 6-MAM                 | < 0.25                              | < 0.2                               |
| Cocaine               | < 0.1                               | < 0.05                              |
| Norketamine           | < 0.2                               | < 0.1                               |
| EDDP                  | < 0.15                              | < 0.1                               |
| Zaleplone             | < 0.2                               | < 0.15                              |
| Zopiclone             | < 0.1                               | < 0.1                               |
| Norbuprenorphine      | < 0.5                               | < 0.4                               |
| Ketamine              | < 0.05                              | < 0.05                              |
| Nitrazepam            | < 0.25                              | < 0.25                              |
| Flunitrazepam         | < 0.5                               | < 0.2                               |
| Clonazepam            | < 0.5                               | < 0.5                               |
| α-OH-triazolam        | < 0.5                               | < 0.25                              |
| Oxazepam              | < 0.2                               | < 0.2                               |
| Estazolam             | < 0.2                               | < 0.1                               |
| Temazepam             | < 0.1                               | < 0.1                               |
| Zolpidem              | < 0.1                               | < 0.1                               |
| Alprazolam            | < 0.15                              | < 0.1                               |
| Methadone             | < 0.15                              | < 0.1                               |
| Lorazepam             | < 0.75                              | < 0.5                               |
| α-OH-alprazolam       | < 0.5                               | < 0.25                              |
| 2-OH-ethyl-flurazepam | < 1                                 | < 1                                 |
| Triazolam             | < 0.2                               | < 0.1                               |
| Nordiazepam           | < 0.2                               | < 0.1                               |
| Diazepam              | < 0.2                               | < 0.1                               |
| Midazolam             | < 0.1                               | < 0.1                               |
| Fentanyl              | < 0.15                              | < 0.1                               |
| Flurazepam            | < 0.2                               | < 0.1                               |
| Buprenorphine         | < 0.5                               | < 0.35                              |

## Additional Notes

### Buffer Preparation

1. 100 mM Ammonium acetate buffer at pH5: Dissolve 7.708 g ammonium acetate in H<sub>2</sub>O and make up to 1L. Formic acid (concentration 98–100% Sigma Aldrich) was used to adjust the pH to 5.
2. 2 mM ammonium acetate aq: Weigh 0.15416 g and dissolve in H<sub>2</sub>O. Dilute and make up to 1 L in H<sub>2</sub>O .
3. 2 mM ammonium acetate in MeOH: Weigh 0.15416 g and dissolve in H<sub>2</sub>O. Dilute and make up to 1 L in MeOH.
4. Dichloromethane/isopropanol (95/5, v/v): Take 95 mL of dichloromethane and add 5 mL of isopropanol
5. Reconstitution buffer: Take 80 mL of 2 mM ammonium acetate (aq) and add 20 mL of 2 mM ammonium acetate in MeOH

### Blowdown Stability

Amphetamines, bath salts and ketamines suffer blow down issues when drying in the free base form. To combat this effect we added 100 µL of 50 mM HCl in MeOH to the collection plate/culture tubes to convert to the corresponding HCl salt forms.

**Reconstitution volumes can be reduced in order to reach lower limits of quantitation, if required!**

## Ordering Information

| Part Number  | Description   | Quantity |
|--------------|---|----------|
| 820-0200-P01 | ISOLUTE SLE+ 200 Supported Liquid Extraction Plate    | 1        |
| 820-0140-C   | ISOLUTE SLE+ 1 mL Supported Liquid Extraction Columns | 30       |
| 121-9600     | VacMaster-96 Sample Processing Manifold               | 1        |
| PPM-96       | Biotage PRESSURE+ 96 Positive Pressure Manifold.      | 1        |
| SD-9600-DHS  | SPE Dry sample evaporator                             | 1        |
| C103264      | TurboVap 96   | 1        |

#### EUROPE

Main Office: +46 18 565900  
Toll Free: +800 18 565710  
Fax: +46 18 591922  
Order Tel: +46 18 565710  
Order Fax: +46 18 565705  
order@biotage.com

#### NORTH AMERICA

Main Office: +1 704 654 4900  
Toll Free: +1 800 446 4752  
Fax: +1 704 654 4917  
Order Tel: +1 704 654 4900  
Order Fax: +1 434 296 8217  
ordermailbox@biotage.com

#### JAPAN

Tel: +81 3 5627 3123  
Fax: +81 3 5627 3121  
jp\_order@biotage.com

#### CHINA

Tel: +86 21 2898 6655  
Fax: +86 21 2898 6153  
cn\_order@biotage.com

To locate a distributor,  
please visit our website at  
[www.biotage.com](http://www.biotage.com)