

# Extraction of Barbiturates from Oral Fluid Using ISOLUTE® SLE+ after Collection with the Oral-Eze® Collection Device Prior to GC/MS Analysis

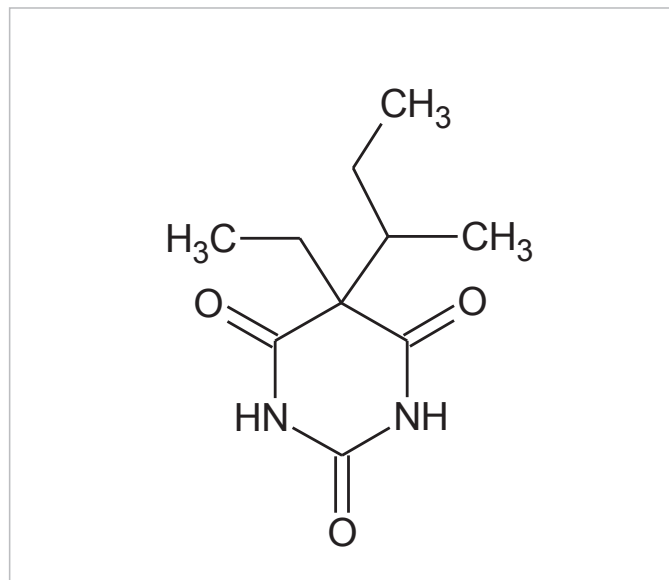


Figure 1. Structure of Butabarbital

## Introduction

This application note describes the extraction of Butalbarbital, Butabarbital, Amobarbital, Pentobarbital, Secobarbital, Hexobarbital and Phenobarbital from oral fluid matrix collected using the Oral-Eze® collection devices, prior to GC/MS analysis.

ISOLUTE® SLE+ Supported Liquid Extraction plates and columns offer an efficient alternative to traditional liquid-liquid extraction (LLE) for bio-analytical sample preparation, providing high analyte recoveries, no emulsion formation, and significantly reduced sample preparation.

This application note describes an effective and efficient ISOLUTE SLE+ protocol optimized for the 400 µL sample capacity column format. The simple sample preparation procedure delivers clean extracts and analyte recoveries greater than 95% with RSDs lower than 9% for all analytes.

## Analytes

Butalbarbital, Butabarbital, Amobarbital, Pentobarbital, Secobarbital, Hexobarbital and Phenobarbital

## Sample Preparation Procedure

- Sample pre-treatment:** Following collection, add 4% ammonium hydroxide (aq) (10 µL) to each collection device (see notes for buffer preparation instructions).
- Format:** ISOLUTE® SLE+ 400 µL Sample Volume Columns, part number 820-0055-B
- Sample loading:** Load 300 µL of the pre-treated oral fluid onto the column and apply a pulse of vacuum or positive pressure (3–5 seconds) to initiate flow. Allow the sample to absorb for 5 minutes.
- Analyte Extraction:** Apply dichloromethane/isopropanol (95/5, v/v, 1 mL) and allow to flow under gravity for 5 minutes. Apply a further aliquot of DCM/IPA (95/5, v/v, 1 mL) and allow to flow for another 5 minutes under gravity. Apply vacuum or positive pressure (5–10 seconds) to complete elution.
- Post Elution & Reconstitution:** Dry the extract in a stream of air or nitrogen using a SPE Dry (40 °C, 20 to 40 L/min) or TurboVap® (1.0 bar at 40 °C for 40 mins).
- Upon dryness, reconstitute with 80 µL ethyl acetate and 20 µL TMAH (trimethylanilinium hydroxide 0.2M) and vortex for 20 seconds. Transfer to a high recovery glass vial.

## GC Conditions

<b>Instrument:</b>	Agilent 7890A with QuickSwap
<b>Column:</b>	Phenomenex Zebron ZB-Semivolatiles, 30 m x 0.25 mm ID x 0.25 µm
<b>Carrier</b>	Helium 1.2 mL/min (constant flow)
<b>Inlet:</b>	150 °C, Splitless, purge flow: 50 mL/min at 1.0 min
<b>Injection:</b>	1 µL
<b>Wash solvents:</b>	Ethyl acetate
<b>Oven:</b>	Initial temperature 120 °C, hold for 1 minute Ramp 12 °C/min to 192 °C Ramp 120 °C/min to 330 °C, hold for 0.85 minutes
<b>Post run:</b>	Backflush for 2.4 minutes (3 void volumes)
<b>Transfer Line:</b>	280 °C

## MS Conditions

<b>Instrument:</b>	Agilent 5975C
<b>Source:</b>	230 °C
<b>Quadrupole:</b>	150 °C
<b>MSD mode:</b>	SIM


## SIM Parameters

**Table 1.** Ions acquired in the Selected Ion Monitoring (SIM) mode

SIM Group	Analyte	Target (Quant) Ion	1 <sup>st</sup> Qual Ion	2 <sup>nd</sup> Qual Ion
1	Butalbarbital	196	195	181
1	Butabarbital	169	184	211
2	Amobarbital	169	184	225
2	Pentobarbital	169	184	225
3	Secobarbital	196	195	181
4	Hexobarbital	235	81	169
4	Phenobarbital	232	146	175

## Results

The optimized ISOLUTE® SLE+ protocol demonstrated analyte recoveries ranging from 95–101% as shown in **Figure 2**. RSDs were below 9% for all analytes.

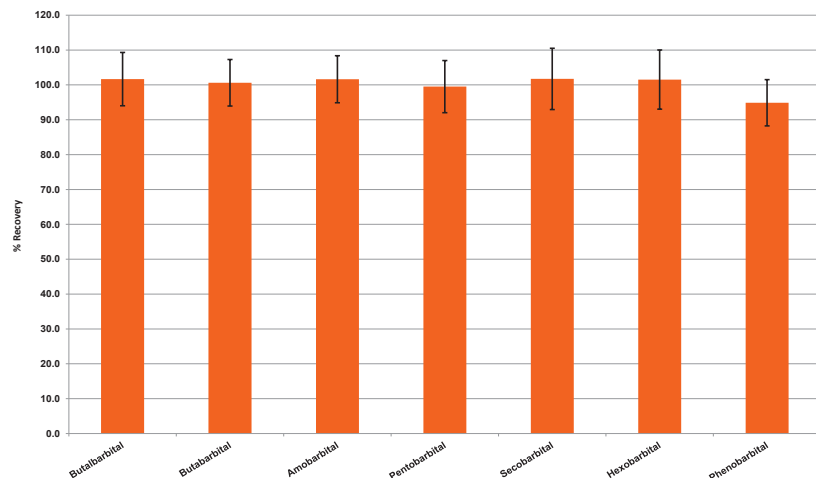

**Figure 2.** Typical extraction % recoveries (n=7) using the ISOLUTE® SLE+ protocol, at a concentration of 500 ng/mL of oral fluid.

## Additional Information

If a non-chlorinated elution solvent is preferred, MTBE (methyl-tert-butyl-ether) is a suitable substitute solvent.

4% ammonium hydroxide is prepared from concentrated stock (28–30%) by adding 200 µL to 5 mL HPLC grade water.

Pre-treating to basic conditions accommodates extraction of multiple basic drugs of abuse alongside the barbiturates class of drugs, in a single assay, if required.

## Ordering Information

Part Number	Description	Quantity
820-0055-B	ISOLUTE® SLE+ 400 µL Supported Liquid Extraction Columns	50
PPM-48	Biotage® PRESSURE+ 48 Positive Pressure Manifold 4	1
SD-9600-DHS-EU	Biotage® SPE Dry Sample Concentrator System 220/240 V	1
SD-9600-DHS-NA	Biotage® SPE Dry Sample Concentrator System 100/120 V	1
C103198	TurboVap® LV, 100/120V	1
C103199	TurboVap® LV, 220/240V	1

For the latest application notes visit [www.biotage.com](http://www.biotage.com)

### EUROPE

Main Office: +46 18 565900  
 Toll Free: +800 18 565710  
 Fax: +46 18 591922  
 Order Tel: +46 18 565710  
 Order Fax: +46 18 565705  
[order@biotage.com](mailto:order@biotage.com)  
 Support Tel: +46 18 56 59 11  
 Support Fax: +46 18 56 57 11  
[eu-1-pointsupport@biotage.com](mailto:eu-1-pointsupport@biotage.com)

### NORTH & LATIN AMERICA

Main Office: +1 704 654 4900  
 Toll Free: +1 800 446 4752  
 Fax: +1 704 654 4917  
 Order Tel: +1 704 654 4900  
 Order Fax: +1 434 296 8217  
[ordermailbox@biotage.com](mailto:ordermailbox@biotage.com)  
 Support Tel: +1 800 446 4752  
 Outside US: +1 704 654 4900  
[us-1-pointsupport@biotage.com](mailto:us-1-pointsupport@biotage.com)

### JAPAN

Tel: +81 3 5627 3123  
 Fax: +81 3 5627 3121  
[jp\\_order@biotage.com](mailto:jp_order@biotage.com)  
[jp-1-pointsupport@biotage.com](mailto:jp-1-pointsupport@biotage.com)

### CHINA

Tel: +86 21 2898 6655  
 Fax: +86 21 2898 6153  
[cn\\_order@biotage.com](mailto:cn_order@biotage.com)  
[cn-1-pointsupport@biotage.com](mailto:cn-1-pointsupport@biotage.com)

To locate a distributor, please visit our website at [www.biotage.com](http://www.biotage.com)

### Part Number: AN824

© 2014 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. Product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe.

For more information visit [www.biotage.com](http://www.biotage.com).