# Biotage® Lysera

en Unpacking Instructions

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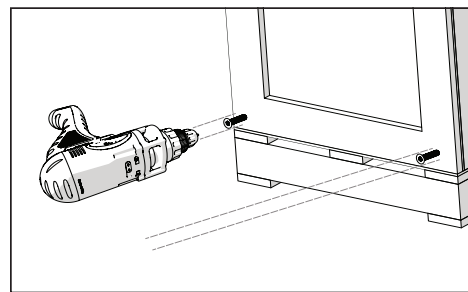
Part Number: UI368.V.1.


© 2017 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. A list of all trademarks owned by Biotage AB is available at [www.biotage.com/legal](http://www.biotage.com/legal). Other product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe.


en

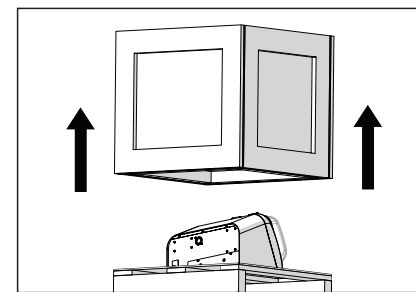
1. Using a screwdriver or power drill, remove the 16 screws around the bottom edge of the crate.

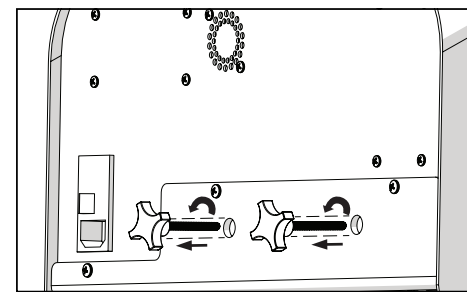
2. Lift the lid from the crate.

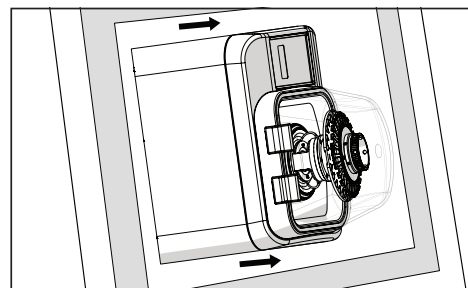
3. Unscrew the two external knob screws from the back of the unit to release from the support plate.  
**WARNING - DO NOT REINSTALL KNOB SCREWS**

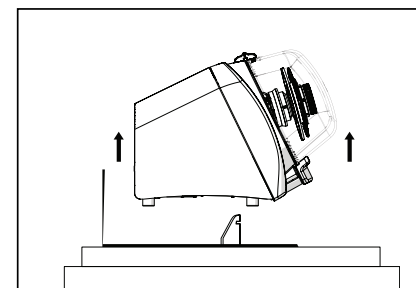4. Slide the Biotage® Lysera away from the knob screw holes.

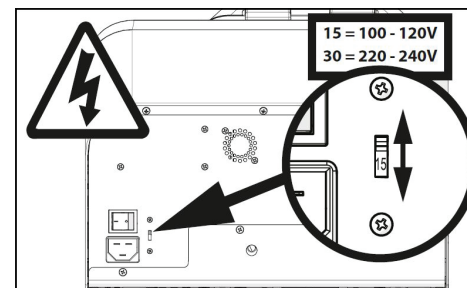
5. Lift the unit up and off of the crate bottom.  
**DO NOT LIFT THE UNIT BY HOLDING THE LID.**

6. **WARNING:** Remove the warning label and make sure the voltage is set correctly for your location before connecting the power cord. If correct voltage is not displayed, use a small screwdriver to slide the red switch up or down to select the correct power supply. Select “15” for 110-120V or “30” for 225-240V power supplies.

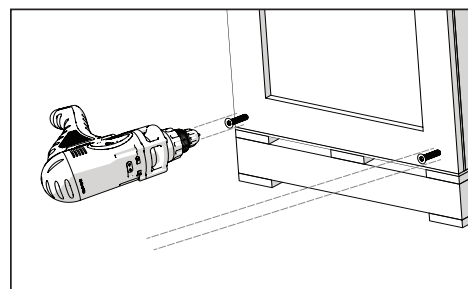NOTE: Do not discard the crate & packaging foam.  
This packaging must be used in the event the Biotage® Lysera needs to be returned for any reason.  
SHIPPING THE BIOTAGE® LYSERA IN ANY OTHER PACKAGING WILL VOID ALL WARRANTIES.

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