

ASE Compatible TurboVap® II

Designed to Accommodate the Dionex ASE™ 40 mL Stemmed Tubes

How Does It Work?

The TurboVap II evaporation system concentrates organic solvent extracts using steam and nitrogen. An optical sensor allows for unattended operation, by stopping the concentration at the 1 mL endpoint. This set-up is ideal for environmental analysis. The patented vortex shearing action provides rinsing as the sample concentrates to maximize recoveries. The six position standard internal rack in the TurboVap II can be replaced with a six position rack to accommodate the new 40 mL stemmed and graduated Dionex ASE vials.

How Will It Benefit My Lab?

With the Dionex ASE and the New ASE Compatible TurboVap II, extractions that previously involved much glassware and hours of manual labor can now be completed with a few minutes of manual labor and one piece of glassware. Productivity is therefore increased and consistency between samples is established. Some of the other advantages include safety, minimized solvent use, minimized sample handling, and no hood space required for either the ASE or the TurboVap.

Can I Convert My Current Turbovap II To be Dionex ASE compatible?

Yes. This new TurboVap configuration is available in two forms. It is offered as a complete TurboVap II already configured for the new stemmed Dionex ASE tube, or as an upgrade kit for current TurboVap II users to convert their current configuration.

Ordering Information

Part Number	Description
C103196	ASE Compatible TurboVap II 110v
C103197	ASE Compatible TurbVap II 220v
C64676	TurboVap II ASE Upgrade Kit

Note: The New Dionex ASE 40 mL graduated tube requires an adaptation to the ASE carousel. This tube is available through Dionex Corporation. Please see www.dionex.com for contact information.

Contact your local Biotage representative for further information. See www.biotage.com for contact information.

NORTH AMERICA

Main Office: +1 704 654 4900
Toll Free: +1 800 446 4752
Fax: +1 704 654 4917
Order Tel: +1 704 654 4900
press (4) at the auto attendant
Order Fax: +1 434 296 8217
ordermailbox@biotage.com

EUROPE

Main Office: +46 18 56 5900
Fax: +46 18 59 1922
Order Tel: +46 18 56 57 10
Order Fax: +46 18 56 57 05
order@biotage.com

JAPAN

Tel: +81 422 28 1233
Fax: +81 422 28 1236
jp_order@biotage.com

To locate a distributor please
visit our web site at
www.biotage.com.


Biotage
www.biotage.com

© 2010 Biotage. All rights reserved. All brand and product names are trademarks or registered trademarks of their respective companies. The information contained in this document is subject to change without notice.