

Operating Instructions

Type 90 and Type 3000 Switching Valves

1.0 DESCRIPTION

Type 90 and Type 3000 valves are manual switching valves containing six or seven ports for a wide range of applications. Type 90 ports accept 1/16" fittings. Type 3000 ports accept 1/8" fittings and are for preparative scale applications. All Type 90 valves, Models 3000, and 3030 are PEEK versions, while Models 3000-038 and 3030-038 are stainless steel versions. Model 9010 contains a position sensing switch.

Model 9010 can be used as a manual sample injector with two-positions. A sample loop attached at Port 1 and Port 4 comes standard. When the sample loop is repositioned to Port 1 and Port 5, Model 9010 can be used in a 4-Way configuration. Removing the sample loop completely makes Model 9010 into a switching valve.

Model 3000 or 3000-038 can also be used as a 4-Way valve (two-position) by connecting a loop to Port 1 and Port 5. Models 9030, 3030, and 3030-038 are Double 3-Way valves (two-position) and Model 9060 is a Six-Position valve.

Figure 1 shows the flow diagrams of Type 90 and Type 3000 switching valves. The circles represent the ports in the valve stator. The dark grooves represent the connecting passages in the rotor seal. See Figure 3 for the flow diagram of Model 9010 as an injector.

2.0 SUPPLIED WITH THE VALVE

Supplied in a separate bag are RheFlex® Fittings sets for all ports and the following items. Model 9010 also includes a 20 μ L sample loop in a separate bag.

- Hex Keys
- Mounting Screws
- Luer Tube (Model 9010 only)

3.0 SPECIFICATIONS

- Maximum Operating Pressure: Models 3000 and 3030 - 28 MPa (276 bar, 4000 psi); All other models - 34 MPa (345 bar, 5000 psi)
- Maximum Operating Temperature: 50°C
- Flow Passage Diameters: Type 90 - 0.3 mm (0.013") and 0.5 mm (0.018"); Type 3000 - 1.0 mm (0.040")
- Wetted Surfaces: Models 9010, 9030, 9060 - PEEK, ceramic, and an inert polymer; Models 3000-038, 3030-038 - stainless steel and PEEK; Models 3000, 3030 - PEEK

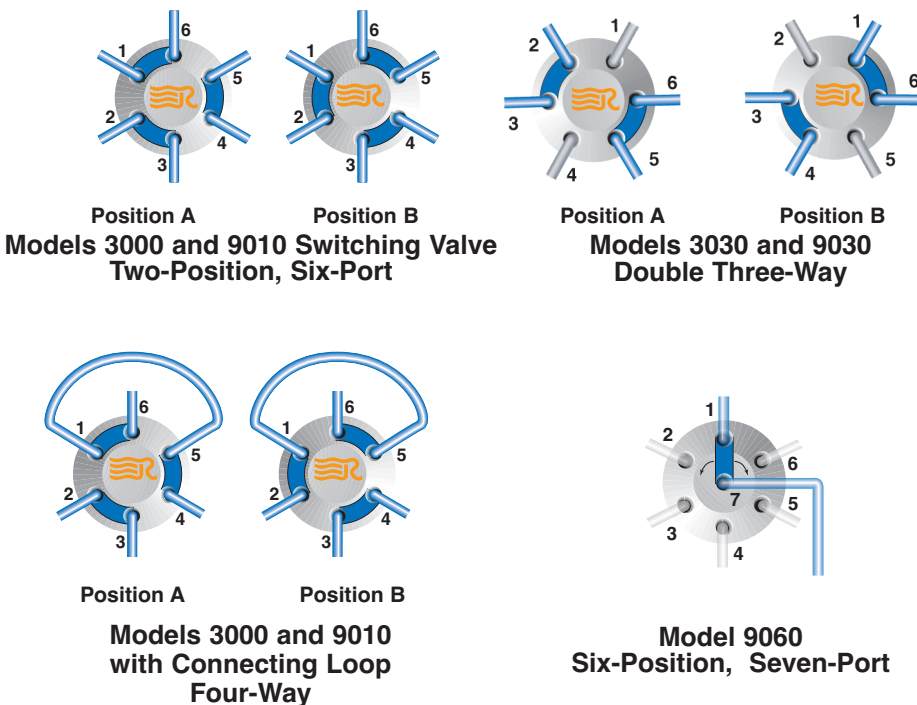


Fig. 1. Flow diagrams of IDEX Health & Science switching valves. The stainless steel versions have the same flow diagrams.

4.0 IMPORTANT SAFETY NOTICES

4.1 Warning: (Using 9010 as an injector): When using sample loops larger than 100 μ L, shield yourself from mobile phase coming out of the needle port when the valve is turned from INJECT to LOAD. Example: 1 mL loop ejects 20 μ L upon decompression from 19 MPa (200 bar, 2898 psi).

4.2 Warning: Burst pressure of PEEK tubing depends on dimensions, solvents, temperature, and duration of exposure. A pinhole rupture in PEEK tubing can emit a high velocity jet of fluid that can penetrate skin and eye tissue.

4.3 Caution: Do not mount the valve with the ports facing up. Leakage due to a damaged rotor seal or loose fittings can cause the bearings to corrode.

4.4 Caution: Rinse the valve thoroughly after using buffer solutions to prevent salt crystals from forming, which can cause damage to the rotor seal and stator face assembly.

4.5 Caution: When using PEEK valves, use only RheFlex® PEEK fittings sets in the stator ports. Metal ferrules can cause irreparable damage to the PEEK stator.

5.0 INSTALLATION

All switching valves can be panel mounted. Model 9010, as a sample injector, connects to different system components. See **Caution 4.5**.

a) Connect the luer tube to Port 5 and the vent line to Port 6. Place the injection port at the same horizontal level as Vent Line 6 to avoid siphoning.

b) Connect the pump to Port 2 and the column to Port 3. Leave the column disconnected from the valve during initial flushing.

6.0 OPERATION

6.1 INJECTIONS (MODEL 9010 ONLY)

Before connecting the column to the injector, flush the injector with mobile phase in both the LOAD and INJECT positions. After flushing the injector, turn to LOAD, and connect the column.

With the accessory injection port (P/N 9012 or 9013) or the luer tube supplied, you can load the injector by two methods: Complete-Filling and Suction Loading.

6.1.1 COMPLETE LOOP FILLING

In complete-filling, the volume of sample injected is set by the volume of the loop (this includes the valve passages). This method produces the highest precision.

Overfill the loop with at least two to five loop volumes of sample. Six to ten loop volumes will provide even better precision. An excess of sample is needed because mobile phase near the wall of the loop is displaced slowly due to the laminar flow effect shown in Figure 2.

To completely fill the loop (see Figure 3):

- See **Warning 4.1** and turn to **LOAD**.
- Insert the syringe into the injection port.
- Load the sample.
- Leave the syringe in position and turn to **INJECT**.

6.1.2 SUCTION LOADING

The preceding method exposes the sample to the metal needle of the loading syringe. Metal can be completely avoided by using a syringe to draw sample into the loop. The steps are as follows (see Figure 4):

- In **LOAD**, dip the tube attached to Port 6 into the sample.
- Insert an empty syringe into the injection port and draw sample into the loop.
- Leave the syringe in position and turn to **INJECT**.

The metal needle of the syringe will contact the sample if an excess of sample is drawn into the syringe, but this excess sample is external to the sample loop and will not be injected. The syringe can be used many times before it needs to be emptied.

To load the loop with all of the available sample, the loop should be at least four times the volume of sample loaded. The loop is first filled with mobile phase via the dip tube, then the whole sample is drawn into the dip tube and loop, followed by more mobile phase. The sample is now sandwiched between two zones of mobile phase in the loop.

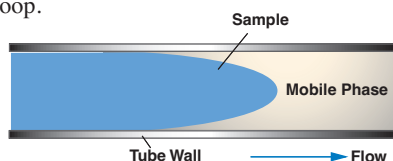


Fig. 2. Laminar flow effect.

6.2 MODEL 9010 AS A 4-WAY

Model 9010 can serve as a switching valve and a 4-Way valve by simply removing or changing the position of the external sample loop (see Figure 1). For 4-Way operation, reposition the 20 μ L sample loop to Ports 1 and 5.

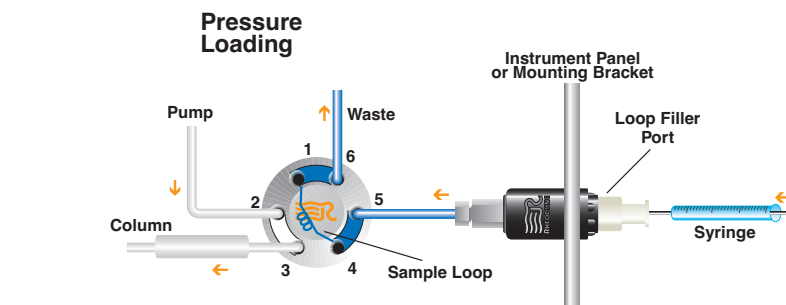


Fig. 3. Using a luer tube to fill the loop by pressure loading.

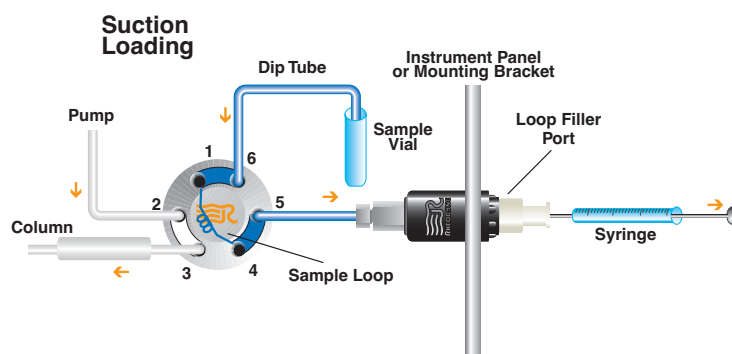


Fig. 4. Using a luer tube to fill the loop by suction loading. In suction loading sample is sucked from a vial into the loop using a dip tube.

6.3 SUGGESTED APPLICATIONS FOR SWITCHING VALVES

Flush the valve with mobile phase before connecting the valve to system components.

Suggested applications for switching valve applications are shown in Figures 5, 6, and 7.

7.0 ADJUSTING FOR LEAKAGE OR HIGHER PRESSURE OPERATION

A single pressure adjusting nut at the handle end of the valve is used for pressure adjustment. If you need operation up to a higher pressure or if there is a leak between the stator and stator ring, remove the handle set screws. This gives access to the pressure adjusting nut. Using a wrench, tighten the adjusting nut about 1/20th turn. Use the red dot and punch mark as guides.

If there is still leakage at this new setting, repeat the process. Finish by retightening the handle set screw into the hole on the shaft. Replace the rotor seal and stator face assembly if the leak continues.

Note: When the valve is not panel mounted, the pressure adjusting nut can be hard to turn. In this case, loosen the three stator screws a 1/4 turn before adjusting the nut. Retighten the stator screws before testing for leaks.

Note (Model 9010 Injector only): Vent Line 6 should be the same horizontal level as the luer tube to avoid siphoning. A siphoning leak will stop when the vent line and sample loop are empty. A leak due to a damaged rotor seal will continue.

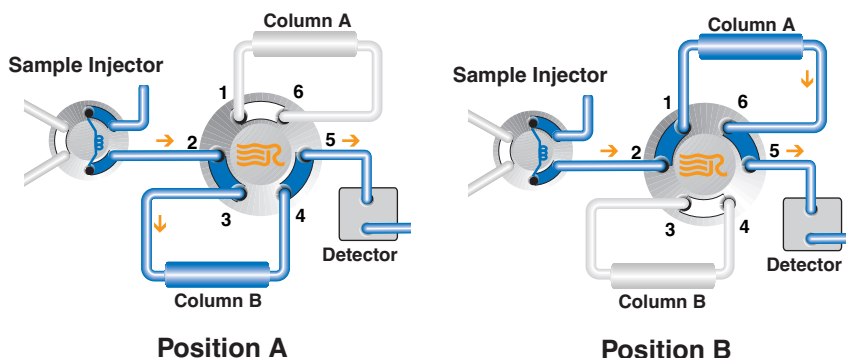


Fig. 5. Two column selection using Model 3000, 3000-038, or 9010 (without the loop).

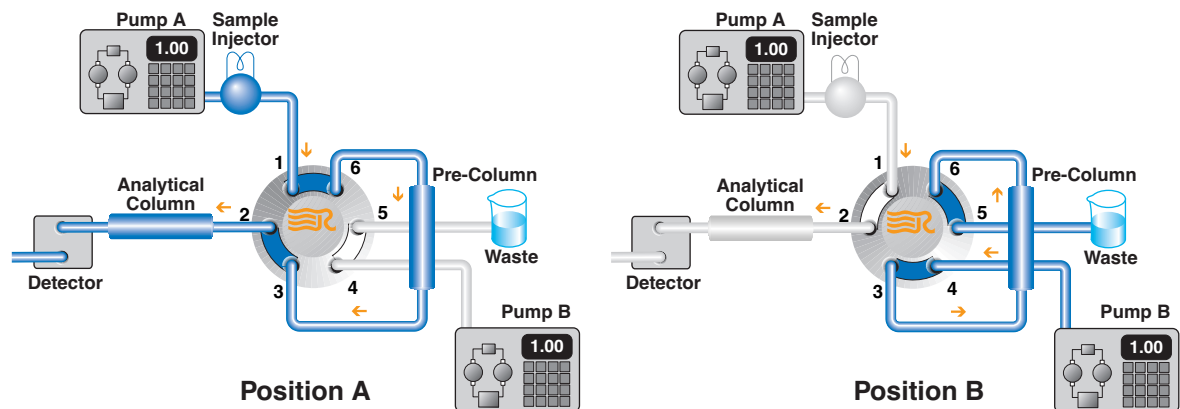


Fig. 6. Stripping method of sample clean-up using Model 3000, 3000-038, or 9010 (without the loop).

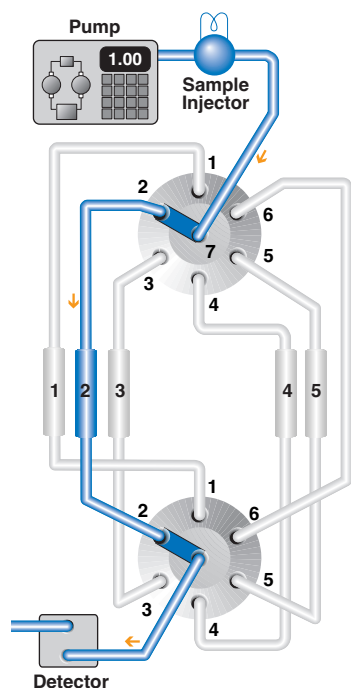


Fig. 7. Five column selection using two Model 9060 valves. The tube connecting Port 6 in the valves is used as a flush-out line.

8.0 MAINTENANCE

The only parts that may need eventual replacement are the rotor seal and stator face assembly. Abrasive particles in the sample can damage the rotor seal and stator face assembly surfaces.

Genuine IDEX Health & Science® parts are easily replaced by the following instructions.

8.1 DISASSEMBLY

To disassemble the valve, refer to Figure 8 and proceed as follows:

- Remove the handle assembly by loosening the handle set screw(s).
- Remove the valve if mounted on a panel.
- Remove the three stator screws.
- Remove the stator support ring (Type 90

valves), stator, stator face assembly, and stator ring from the body.

e) Pull the rotor seal off the pins.

f) Leave the isolation seal and bearing ring in place. On Model 9010, remove the isolation seal.

8.2 REASSEMBLY

To reassemble the valve, refer to Figures 8 and 9 and proceed as follows:

a) On Model 9010, put the new isolation seal on the shaft assembly with the open side facing the handle.

b) Orient the new rotor seal as shown in Figure 9. For six-position valves, orient the rotor seal on the rotor relative to the arrow on the shaft. The rotor seal grooves face the stator.

c) Replace the stator ring so that the pin in the stop ring enters the mating hole in the stator ring. For Type 90 valves, confirm that the position sensing switch falls into the mating hole in the stator ring.

d) On Type 90 valves, place the stator support ring onto the stator. Be sure the stator support ring's bumpy inner contour fits into the mating contour of the stator.


e) Place the pins in the new stator face assembly into the mating holes in the stator. The three stator face assembly pins in the Type 3000 valves fit into the mating holes in the stator only one way.

f) Replace the stator and stator face assembly on the valve so that the pin in the stator ring enters the mating hole in the stator.

g) Replace the three stator screws. Tighten each screw a 1/2 turn past fingertight.

h) If panel mounted, remount the valve onto the panel.

i) Replace the handle assembly and tighten

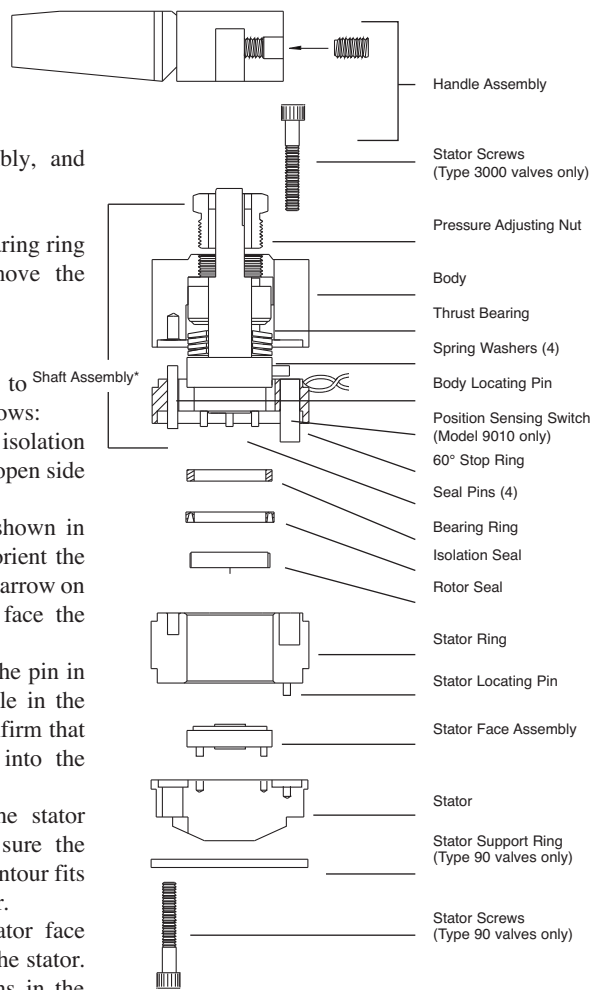

* Shaft assembly includes shaft, rotor, and pins.

Fig. 8. Exploded view of IDEX Health & Science switching valves.

the two handle set screws into the hole on the shaft.

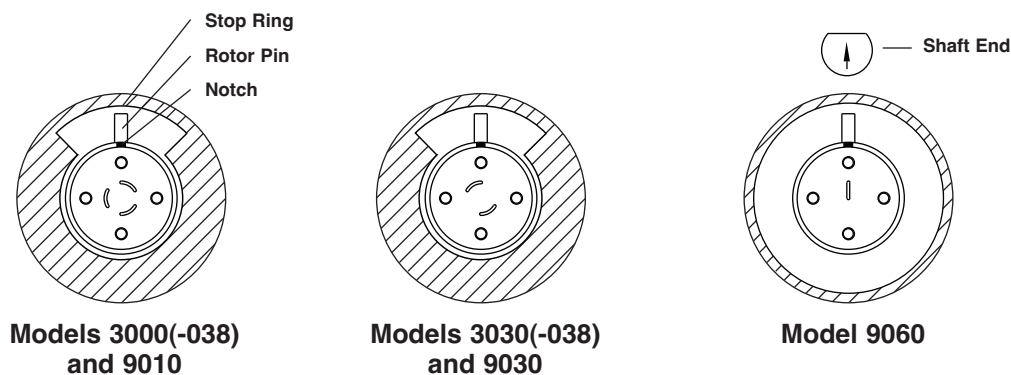


Fig. 9. Correct rotor seal orientation as viewed from the stator.

8.3 POSITION SENSING SWITCH IN MODEL 9010

The switch is a magnetic reed switch actuated by a magnet sealed inside the shaft. The switch is rated for 100 V at 200 mA.

To replace or remove the switch:

- Remove the three stator screws.
- Remove the stator support ring, stator, stator ring, and stop ring.
- Pull the switch out of the stop ring.
- Replace with new switch, or leave the hole empty if the switch is not needed.
- Follow the steps in Section 8.2 to reassemble.

8.4 DETENT MECHANISM (SIX-POSITION VALVES ONLY)

If mounting the six-position valve on an actuator, the detent mechanism must be disabled:

- Remove the stator screws, stator support ring, stator, stator ring, and stop ring.
- Loosen the handle set screw(s) on the knob and remove the handle assembly.
- Push the shaft through the body.
- Remove the two springs and two balls from the stop ring.
- Replace the shaft back into the body (shaft fits only one way).
- Follow the steps in Section 8.2 for reassembly.

9.0 OPERATING SUGGESTIONS AND TROUBLESHOOTING

9.1 LEAKAGE

If you see liquid between the stator and stator ring, tighten the pressure adjusting nut as explained in Section 7.0. Replace the rotor seal and stator face assembly if the leak continues.

9.2 USE OF AQUEOUS BUFFERS OR SALT SOLUTIONS

To prevent the formation of salt crystals in the valve, flush out the passages and vent line(s) with water after using salt solutions.

9.3 ACCURACY OF SAMPLE LOOPS

Sample loop sizes are not actual values. The actual volume can differ by $\pm 10\%$ for a 20 μL loop and $\pm 5\%$ for a 10 mL loop. There is a greater difference for smaller loops.

10.0 RECOMMENDED SPARE PARTS

IDEX Health & Science offers a RheBuild® Kit, including all the necessary parts, tools, and instructions for Model 9010, and offers genuine IDEX Health & Science® replacement parts to maintain the quality performance of your valve.

9010-999 RheBuild Kit for Model 9010

7030-015 Rotor Seal for Model 9030

7060-074 Rotor Seal for Model 9060

3710-008 Rotor Seal for Models 3000 and 3000-038

3030-005 Rotor Seal for Models 3030 and 3030-038

8125-094 Stator Face Assembly for Model 9030

9060-015 Stator Face Assembly for Model 9060

11.0 WARRANTY

All IDEX Health & Science products are warranted against defects in materials and workmanship for a period of one year following the date of shipment by IDEX Health & Science. IDEX Health & Science will repair or replace any IDEX Health & Science product that fails during the warranty period due to a defect in materials or workmanship at no charge to the customer. The product must be returned to IDEX Health & Science's factory in original packaging or equivalent, transportation prepaid. Damage occurring in transit is not covered by the warranty. This limited warranty is IDEX Health & Science's sole warranty of its products, and all other warranties of merchantability or fitness for any particular purpose are hereby disclaimed. Under no circumstances will IDEX Health & Science be liable for any consequential or incidental damages attributable to a claimed failure of a IDEX Health & Science product, even if IDEX Health & Science has been placed on notice of possibility of such damages.